## D

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#### Batterie

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### Operating Instructions

#### Battery

An 9V block battery is included; this should be placed into the battery compartment in the manner indicated on the base plate.

#### Positioning of the Scales

Please ensure that the scales are in a horizontal position.

### Weighing (🔊)

After switching on the scales with the **🔊**-button, all segments are shown on the display. Please wait until the zero appears. Then place the weight on the scales and read off the weight shown.

#### Net Weighing (tare)

Place an empty container (or the first weight) on the scales and press the „tare“-button until the zero appears. Fill the container (or place the second weight on the scales). Only the additional weight is indicated in the display.

### Switch-off (🔊)

Press the **🔊**-button. Automatic switch-off after 2 minutes whether weight is on the scales or not, if no change of weight occurs in this time.

**Changing the weighing units: Kg, Lb (mode)**
This scales can show the weight in either kg or lbs. Press the „mode“-button until the required weighing unit appears.

### Counting (🔊🔊🔊)

When the „zero“ appears place the reference weight on the scales. Then press the **🔊🔊🔊**-button and hold until „100“ appears in the display. Press the „mode“-button and choose the reference quantity (25; 50; 75 or 100) and the press the **🔊🔊🔊**-button once again.

#### User Calibration

If needed the scales can be recalibrated. Ensure that the scales is turned off and press the “mode” and „tare“-buttons and keep them pressed while the scales is turned on. **🔊**. As soon as a numeric code appears, let go the “mode“-button first and then the “tare“-button. Press the „mode“-button until „ZER0“ appears and then a numeric code. Then place a **120 kg**-calibration weight on the scales. Confirm with the „mode“-button the calibration weight. First „120“ appears in the display and then „PASS“ and finally the normal weighing display. The scales is now calibrated . If „FAIL“ instead of „PASS“ appears then the calibration is not correct and the display disappears.The calibration routine has to be repeated. Important: The scales should not experience any movement or draughts!

**Explanation of the special symbols**
1. Switch-On

After pressing the **🔊**-button all symbols appear. One can check if all segments are shown correctly. The “zero” which then appears shows that the scales is ready for weighing.
2. Negative Weight Display
Press the „tare“-button again.
3. Overload
If the weight on the scales is heavier than the max. weight then „O.Ld“ appears in the display.
4. Power Supply
If „LO“ appears, this means that the battery is used up and has to be replaced.

This device corresponds to the requirements stipulated in 90/384/EEC. Note: Extreme electromagnetic influences e.g. radio unit in the immediate vicinity may affect the displayed values. Once the interference has stopped, the product can once again be used normally. This scales is not legal for trade.

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#### Garantee

This product has been manufactured with the utmost care and by applying the strictest quality guidelines. The material used and the manufacturing process are subject to constant testing.

The guarantee is valid for three years and covers the repair of defects in the production process or in the materials used. The period of guarantee commences with the day of sale to the first user. The payment slip or the purchase bill are proof. Damage attributable to improper handling, mechanical damage or natural wear and tear as well as damage to due the batteries (if supplied) are excluded from the guarantee. In cases of manipulation by parties other than the supplier the guarantee lapses. No liability is accepted for consequential damages.

Complaints in respect of recognisable defects must be reported within fourteen days after the receipt of the goods. Hidden defects should be notified immediately on discovery. Goods which have been justifiably rejected will be repaired. In case we let a reasonable period of grace pass without eliminating the defect the buyer is entitled to withdraw from the contract.

In case of guarantee claims please forward the device with the receipt to the point of sale. Please, provide your name, address and a description of the defect.

In the case of defects occurring outside the guarantee period a repair or replacement subject to payment of the costs is possible. In such cases, please contact again the point of sale.

Nothing in the above conditions excludes the application of provisions of the EU.

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#### Precision:

This device corresponds to requirements stipulated in 90/384/EEC and 89/336/EEC. Every scale has been carefully calibrated and controlled during the production process. The tolerance is " 0,5% " 1 digit ( at Temperatures between +5° and +35° C) Incorrect display values due to damage attributable to improper handling, mechanical damage or malfunction is exempt from liability. Damages due to faults are also exempt from the guarantee. No liability is accepted for consequential damages or losses by the buyer or the user.

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JAKOB MAUL GmbH